

# Installing the FilmArray® BioThreat v2.5 Software Pouch RUO Module

## Introduction

---

In order to use the BioThreat v2.5 pouches with the FilmArray system, the BioThreat v2.5 Software Pouch Module must be installed. **Please note the Software Pouch Module must be installed on a FilmArray laptop running FilmArray Software v1.4.**


Please contact BioFire Customer Support if you have questions about updating to FilmArray Software v1.4.

## Installing the Software Pouch Module

---


Using the included software disk complete the following steps to install the new pouch module on each FilmArray laptop.

1. Insert the BioThreat v2.5 Software Disk or USB Memory Stick into the appropriate drive.
2. Open the FilmArray Instrument Control software from the desktop icon (if not already open).

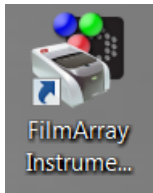


ADVISORY  
::: NOTICE

3. Navigate to the 'FilmArray' menu in the top left of the software. Click 'Tools'. Click 'Pouch Management'.

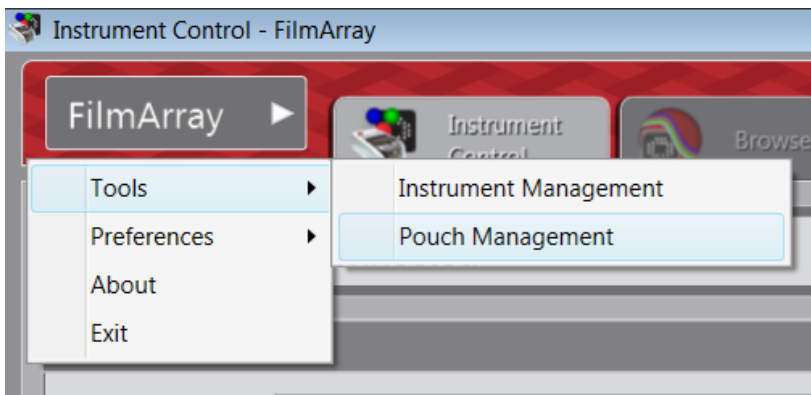

4. Click 'Install New Pouch'.

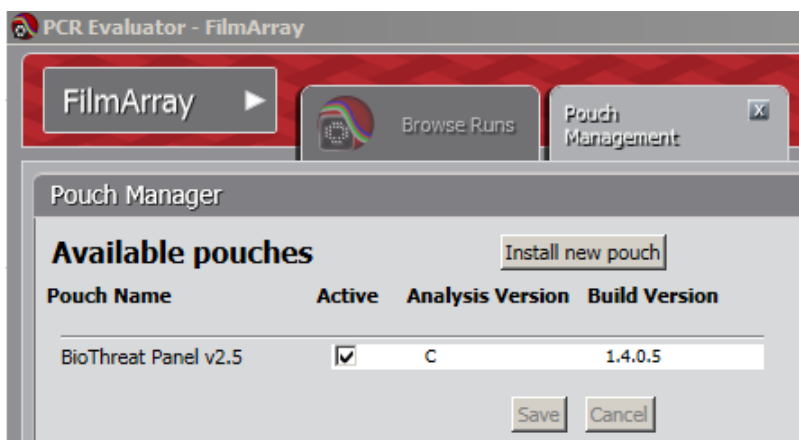


**NOTE:** The list of Available pouches on your system may differ than the picture above.


5. Navigate to the CD / DVD / USB drive in the Open File dialogue.
6. Click the BioThreat Panel v2.5 pouch module file named "BioFireDx.FilmArray.Pouches.BioThreatPanel-2.5.dll".

## ADVISORY NOTICE

7. Click 'Open'.

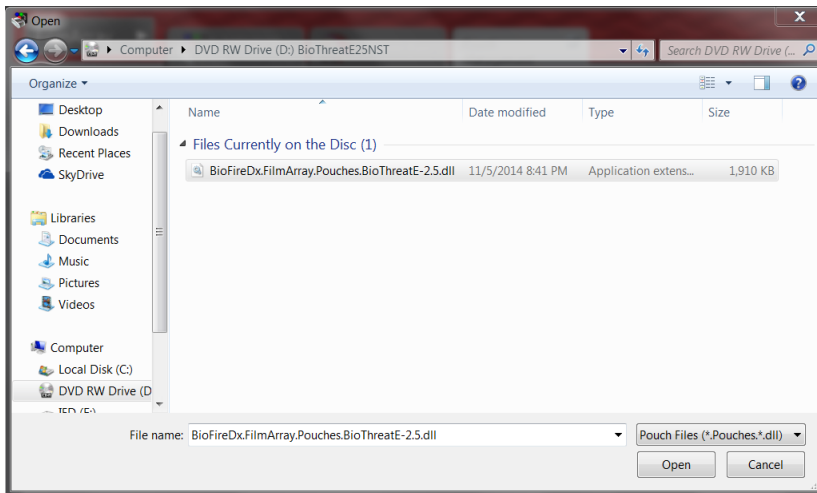

8. The BioThreat v2.5 Pouch Module will display and the 'Active' box will be checked when it has successfully loaded.
9. Verify the BioThreat v2.5 Build version reads 1.4.0.5.

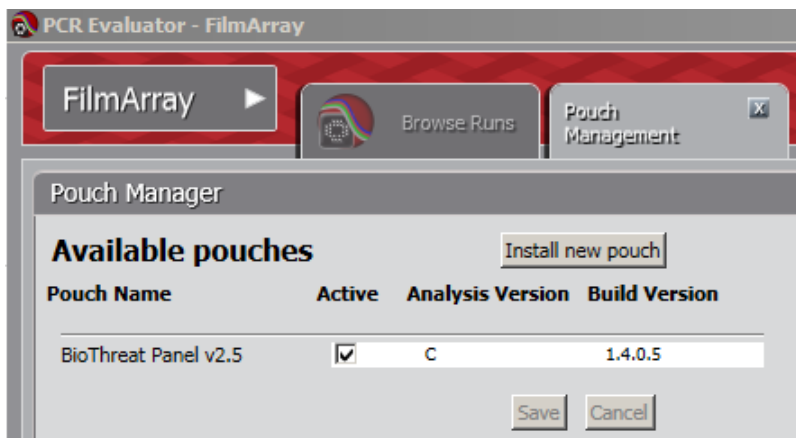

**NOTE:** The list of Available pouches on your system may differ than the picture above, with the exception of BioThreat v2.5.

10. Eject the install disk / USB stick from the laptop.

ADVISORY  
NOTICE

## Technical Support Contact Information

---

BioFire Defense is dedicated to providing the best customer support available. If you have any questions or concerns about this process, please contact the FilmArray Technical Support team for assistance.

### General Information

Email: [support@biofiredefense.com](mailto:support@biofiredefense.com)

Phone: 1-801-262-3592

Fax: 1-801-447-6907

### Kall Symons

Technical Support Specialist

Phone: +1-801-262-3592 extension: 1406

[kall.symons@biofiredefense.com](mailto:kall.symons@biofiredefense.com)



**ADVISORY  
::: NOTICE**